The Product ID and Extra ID tab allows you to set the default product ID/ Extra ID for your standards and samples. A list of Product ID/ Extra ID may be maintained and chosen on their tabs. Type text for a new Product ID/ Extra ID into the box and then click **Add Selection** to add it to the list. Choose an ID to be deleted from the drop down box and click **Remove Selection** to remove it from the list. The ID that is currently selected from the drop-down box is the one that will be automatically applied on reading.

laming Conventions	
Sample Standard Series Proc Product ID list. Type new ID's in Selection'' button to add them to ID and select the "Remove Sele from the list.	to the box and select the "Add the list. Highlight an available
[	<u>.</u>
Add Selection	Remove Selection
Additional Settings Limit the Standard Product Use only existing Product Use only the current Pro- Limit the Sample Product	ct ID's aduct ID selection.
Use only existing Product     Use only existing Product     Use only the current Pro     Set to the Standard's Pro	ct ID's aduct ID selection.
OK.	Cancel Apply

Naming Conventions	
Selection" button to add them	Product ID Extra ID to the box and select the "Add n to the list. Highlight an available selection" button to remove items
Add Selection Additional Settings   Limit the Standard's Ex  Use only existing Ext  Use Extra ID selecte	
<ul> <li>Limit the Sample's Extr</li> <li>Use only existing Ext</li> <li>Use only the current</li> <li>Set to the Standard's</li> </ul>	Extra ID selection.
OK	Cancel Apply