

Why does ColorFlex EZ Tomato or Citrus flash 3 times in EasyMatch QC?

FAQ: "We noticed that CFEZ flashes 3 times per measurement when connected to EasyMatch QC. Do you have any comments on why this might occur?"

When ColorFlex EZ is being used with EasyMatch QC and any of Tomato metrics or Citrus metrics are configured to display in any of Color Data Table/Trend Plot/EZView views, EasyMatch application will prompt for there (3) readings from the ColorFlex EZ and display the average as the measurement.

This is part of CMR implementation as translucent, light trapping orange juice and processed tomato products will benefit from averaging multiple readings to give a more repeatable measurement.